ORTHOLUX II



Instructions




ORTHOLUX II



Instructions

2	Unpacking the microscope												
3	Assembling the microscope												
4	Technical details												
411	FSA binocular tube						14		-				
412	S binocular tube												
42	Objective revolving nosepied	e											
43	Microscope illuminator .												
44	Objectives												
441	Oil immersion objectives												
45	Eyepieces												
5	Operating the microscope												
6	Transmitted-light darkground												
7	Examinations in phase contra	st											
B	Fluorescence												
9	Examinations in incident-ligh	td	ark	gr	out	nd							
10	Microscopic measurement			1									
11	Heating stages												
12	Photomicrography												
	Care and maintenance .												

1 Technical description

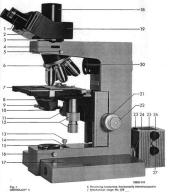

OffTHOLUX* II

1. Widefield eyepieces in the eyepiece tube
2. Moulded grips for the edjustment of the interpupilisty distance

Interpupillary distance
3. Lever for tube change
4. Filter slot
5. Clamping screw for arresting the revolving nosepiece

Swing-out condenser No. 602
 Centring screws on the swing-out condenser No. 602
 Agenture disphrages lever
 Knowled source for the vertical adjustment

of the condenser

12. Co-axial knobs for the adjustment
of the mechanical stage

2 Unpacking the microscope

The following parts are packed in a special container: 1) Microscope stand with changing

guides for the parts to be attached 2) Microscope tube 3) Object stage with changing piece

and condenser carrier 4) Lamp Housing 50

5) Components such as objectives, evenieces, condenser, dust cover,

rial

etc. Transformers and other heavy accessories are packed separately. During unpacking carefully check the equipment against the packing note and make sure that no small components are left among the packing mate-

Touching the lenses of the objectives and evenieces must be avoided if possible. Any fingermarks on glass surfaces must be removed at once with a soft piece of chamois leather or a wellwashed piece of linen. Even slight traces of finger perspiration may attack the surfaces of high-quality optical glasses within a short time

The workroom should be as free as possible of dust, oil vapours and chemical fumes which attack optical and mechanical parts. Nor should there be any great temperature fluctuations or vibra-

tions. The socket for the built-in illuminator should be fused for 13 amp.

13. Knob for the swing-out lens 14. Field disphragm 15. Dust glass with filter support 16. Clamping device for accessories

16. Photo tube with removable dust cap 20. Disphragm tube with filter slot 21. Rotating knob for coarse adjustment

Rotation knot for time ortisatement 24. Locking device for the Lamp Housing 50 25. Lamp Housing 50 25. Knurled knob for the lateral adjustment

27. Knurled knob for the vertical adjustment

3 Assembling the microscope

Release the clamping screw of the object stage (Fig. 229), insert the object stage into the changing guide (228), and lower it until the top of the angle bracket of the stage is flush with the top of the drive casing. Fix the stage with clamping screw (229).



Fig. 2 reserting the object stage

Lower the stage with the coarse adjustment until the objective revolving nosepiece or ULTROPAK® can be inserted fully into the horizontal changing guide (3.30), with the clamping screw (1.5) released. Re-tighten clamping screw (1.5).



Fig. 3 Inserting the revolving nosepiece

Push the lever for the tube change to the rear and insert the tube in the changing bayonet from above. The rotation of the tube is immaterial. Release the locking lever. It must be possible to rotate the tube through 360° freely after insertion. You can clamp the tube by slightly tightening the lever (4.3) towards the front.



Fig. 4 Inserting the FSA binocular tube

Lower the slide changer (5.31) by means of knob (5.11) so that the swing-out condenser can be easily inserted in it as far as it will go. Ensure that the two centring screws face the observer. The rotating knob (5.32) serves for the swinging in and out of the condenser too.

- 11 Knurled knob for raising and lowering the condenser 29 Clamping screw for fixing the object
 - Dovetall changer of the condenser
 Knurled knob for swinging the condenser to
 in and out of the beam

Attaching the Lamp Housing 50 for transmitted-light illumination

Slightly lift the stand or place it at the edge of the table. Turn the Lamp Housing 50 through about 90" to the left, insert it in the bayonet mount, and lock it in position by turning it to the right.



Fig. 5 Inserting the condenser



Fig. 6 Attaching the Lamp Housing 50

Inserting the 12 v 50 W lamp Release the knurled screw (8.33) on the

Lamp Housing 50 and remove the side wall from the housing.
Pull out the 12v 50W lamp (7.35) from the socket. Insert the new lamp with

protective cover in the plug-in socket and remove the protective cover. Insert the side wall so that holder and plugs (7.36) engage. Re-tighten the knurled screw.



Fig. 7 Inserting the 12v 50W lamp in the Lamp Housing 50

Centring the 12v 60 W tungsten halogen lamp in the Lamp Housing 50



Lamp Housing 50

Locking device of the Lamp Housing 50
 Knurled knob for the lateral adjustment of the lamp
 Knurled knob for the vertical adjustment of the lamp
 Knurled knob for the vertical adjustment of the lamp
 Knurled screw for opening the lamp housing



Fig. 9

After each lamp change the tungsten halogen lamp must be re-centred. Fully open the field diaphragm. Place

rully open the field diaphragm. Place the centring disc on the dust glass of the microscope.

a) Focus the mirror image of the lamp filament by adjusting the reflector (8.34).

Form an image of the filament and of its mirror image on the centring disc by rotating ring (8.23).

b) Move the image of the filament and its mirror image into the centre by rotating knob (8.26).

c) Rotate knob (8.27) for the vertical adjustment of the lamp until the image of the filament and its mirror image over-

After insertion of the diffusing disc and removal of the centring disc rotate the focusing knob (8.23) for the lamp condenser, simultaneously observing the image through the eyepiece tube, until the rear focal plane of the objective is evenly illuminated.



Fig. 10 Inserting the 15W lamp in the Lamp Housing 15



Inserting the 6 v 15 W lamp The lamp is changed as follows:

The lamp is changed as follows:

1) Release the arresting screw (10.37) and remove the lamp socket.

2) Remove the lamp (10.38) (push the

lamp into the socket (10.39) and unlock it by turning it to the left). 3) Both pre-focus or non-centred lamps

can be used:

a) insert the pre-focus lamp so that the slot of the lamp lies below the red indicating mark of the lamp socket.

b) insert the non-centred lamp with the two lugs in the guide groove of the socket and lock it in position by pushing it in towards the right.



Insert the Lamp Housing 15 in the bayonet mount, with the arresting screw

for the lamp condenser adjustment pointing about 90° to the left. Lock it by turning it to the right. The arresting knob now points upwards.

Centring the 6v 15 W filament lamp in the Lamp Housing 15

in the Lamp Housing 15
Place the centring disc on the dust glass. Open the field diaphgram. Release the arresting screw and adjust the lamp socket (10.39) until the image of the light point on the centring disc is at its smallers.

Adjust the two centring screws (10.40) until the light points are in the centre of

the centring disc.

After removal of the centring disc adjust the lamp socket until the rear focal plane of the objective is evenly illuminated.

4 Technical details

4 11 FSA binocular tube

The various components such as object stages, tubes, objective revolving nose-pieces, etc. are described in detail in our List ORTHOLUX II. If special descriptions of individual elements are required for correct operation please consult the following paragraphs.



Par supplier to

The FSA binocular tube for the ORTHO-LUX II must be used for all evenieces of 23.2mm diameter. It has a swing-out beam-splitting prism, which either splits the light intensity at a ratio of 80:20 (80% for photomicrography, 20% for visual observation) or directs the entire light (for visual observation) into the eyepiece tubes. The interpupillary distance is set with both hands laterally pulling or pushing the tubes. The optical mechanical-length compensation ensures perfect sharpness for any internunillary distance both in the observation evenieces and in the film plane. Where the interpupillary distance is unknown the tube is adjusted during binocular observation until only one circular and comfortably surveyed field of view appears. Corrections of defective vision must be carried out with the aid of the focusing evelenses of the PERI-PLAN® evenieces.



Fig. 13 5 binocular tube



The tube can be adjusted for the individual interpupillary distance of the observer. This requires an appropriate correction of the tube length, which is carried out on the eyepiece tubes as follows:

Set the interpupillary distance by pushing or pulling with both hands so that the two part images are coincident in the microscope and only one round image is seen. Read the interpupillary distance thus determined off the scale on the front pale of the two bays and trans-

fer it to the two eyepiece tubes.

Example: When an interpupillary distance of 65mm has been determined set the left and right-hand eyepiece tubes at the scale number 65. For defective eyesight one of the eyepiece tubes can be additionally adjusted for sharpness correction.



Fig. 14 Revolving nosepiece

I The objective revolving nosepiece can be be notronatuly exchanged with other of optical elements such as the ULTRO-PAK vertical illiminator. The nosepiece has five numbered threads to receive the objective. Each outfit is accompanied by an optical objective/eyepiece ochart. This indicates, among other data, the eyepiece threads to which the individual objectives are matchetives are matchetives.

43 Microscope illumination

The sharpness and the general character of the image are decisively influenced by the microscope illumination. Besides the choice of correct objectives and reproduction scale careful setting up of the illumination and its adjustment to the specimen are some of the basic conditions for optimum utilization of the

microscope. The following demands must be made of microscope illumination - i.e. the optical path from the light source to the specimen:

1) The illuminator must supply the microscope with a beam cross section required by the objective/evepiece combination in use

2) The illuminator must make it possible to limit this beam cross section at will. Whether and to what extent the first demand is met can be seen from the complete and even illumination of the field of view in the microscope. After removal of the eveniece the exit pupil (generally the rear lens of the objective) must also

appear evenly illuminated at the bottom of the tube

But the maximum beam cross section required by the individual objective/ evepiece combination varies. Maximum diameter is, however, not needed by the highest-power, but by the lowest-power objectives, which cover a larger object field. The condenser must therefore be computed so that it can deliver the beam cross section required by the lowestpower objectives. But this creates a surplus for higher-power objectives which

cannot be fully utilized, and which may even adversely affect image quality through reduction of the contrast, causing reflections, and unnecessary heating of the specimen. It is therefore necessary to be able to reduce the beam cross section to the degree that is applicable and useful for the optical combination used in the microscope. This is achived by two diaphragms suitably arranged in the beam.

The image of one disphragm limits the rays in the specimen. If this diaphragm is closed, it will be visible in the microscope as field diaphragm. It thus offers the possibility of meeting the second demand, in that the light beam serving for the illumination of the field of view can be limited to the desired extent through the opening and closing of the disphragm. The second disphragm serves to meet this demand also for the beam cross section in the east pupil or rear lens of the objective. Its image disphragm. The as a circular boundary of the portion of the rear lens of the objective filled by light after removal of

the eyeplece. For the true Koehler's Illumination realized in the ORTHOLUX II the field diaphragm is situated between the lamp condenser and the deflecting mirror, whereas the aperture diaphragm is built into the condenser. The paragraph "Centring the swing-out condenser" provides information on the use of these

two diaphragms. The achromatic aplanatic swing-out condenser No. 602, N.A. 0.90, is part of the standard outfit of the ORTHO-LUX II. With this aperture the aperture of practically all dry system can be fully illuminated.



Fig. 15 Image of the sporture disphragm in the objective



Condenser bottom part No. 600 with the two condenser tops ACHROMAT 0.90 and APLANAT 1.25

12 Agenture disphragm lever 12 Rotating knob for swinging the condenser to it and out of the beam 41s Condenser top No. 002 41b Condenser top No. 003 Immersion objectives, too, can be used with this condenser; only if the distinction of the most delicate structures is essential will a condenser aperture above 0.90 become necessary. Our

achromatic-aplanatic swing-out condenser No. 603, A 1.25 is available for this purpose. The table below offers information on the properties of the condensers.

Swing-out condensers

	light brightfield Description	Application
600	Bottom part of condenser, A 0.25, with aper- ture diaphragm, condenser lens for low magnification and	For objectives of up to A 0.25. With this condenser even large fields can be illuminated at low magnification.
002	3-lens condenser top Actr. 0.90 forms	
602	Achromatic swing-out condenser Achr. 6.90, of good spherical and chromatic correction or	For more highly corrected objectives, such as fluorite systems, apochromats, plano ob- jectives, and for photomicrography; limited suitability for fluorescence.
003	4-lens condenser top Apl. Oel 1.25 forms	
603	Actromatic-aplanatic swing-out condenser Apl. Oel 1.25, state of correction similar to that of condenser 602 at large apertures.	Mainly for work with highly corrected oil in- mension objectives or for colour photomicro- graphy at large condenser apertures. Un- suitable for fluorescence.

For the investigation of tissue cultures, or very thick specimens etc., condensers of lone intercept distances are used for

Designation	Descr	iption		Working distance	Corrected for		
605	botton	matic condenser No m part 600 and cond- ved Achr. 0.70/L 4	4mm	4mm glass, water Agar-Agar etc.			
606	consis	ohromatized conder sting of bottom part ved 0.60/L 11	ser No. 606, A 0.60, 600 and condenser top 606,	11mm	11mm, of which about 6mm glass etc.		
607	consis	chromatized conder sting of bottom part ved 0.45/1, 20	ser No. 607, A 0.45, 800 and condenser top 007,	20mm	25mm, of which about 6mm glass etc.		

must be seen of the field disphragm.

Lower the swing-out condenser until a sharp image is termed of the field disphragm.

Every microscope objective has a number of data engraved in addition to our firm's emblem, which are important for recognition by the user. The following data are engraved:

170

The distance in mm from the shoulder of the objective to the rim of the tube. This distance is also called mechanical tube length, All LEITZ transmitted-light objectives are corrected for 170mm mechanical tube length. This length cannot be maintained in our tubes with inclined eveniece tube because of the optical elements such as prisms, beam splitters, etc. they contain Nevertheless the objectives are correctly used with such tubes, since a tube lens displaces the image into the new intermediate plane without adverse effect on the image quality. Suitable choice of the tube lens makes it possible to make the image appear at the same reproduction ratio in the intermediate plane



LETZ plano objectives with cross section The PI Apo Cel 100/1.32 objective has 14 elements

0.17

This is the thickness of the coverglass for which most of our transmitted-light to objectives are computed, instead of the figure 0.17 a dash may be engraved. This means that with these objectives speciemen with or without coverglass (e.g., smear preparations) can be observed. The table shows which objectives may be used with or without coverglass (column owners accurations).

Below the details about tube length and coverglass correction the following data will be found in abbreviated form: the reproduction scale (size ratio = intermediate image: object e.g. 10:1) and

the numerical aperture of the objective. here 10303. For reproduction scale the word magnification is used in the table (this denotes objective magnification). In front of the reproduction scale the state of correction with Placetie spatems, Asochorenatis, Neman Place objective or contents, and the placeties of correction with Placeties spatems, Asochorenatis, Objectives for placeties and the placeties of the additional designation Placeties and setting alternative or the additional designation Placeties and the content strength of the placeties that can be used only with an immersion medial internetion of the placeties of the placeties and the placeties of th (PI Apo **Oel** 100/1.32). In addition all immersion objectives display a black ring on their mounts for instant recognition. For the protection of the specimen at the front lens of the objective all the higher-power systems have a springloaded front lens mount.

Objectives for transmitted-light investigation in brightfield or darkground

Designation	of objectives Magnification/aperture	Cover- glass correc- tion 1)	Type of eyepiece	Free work- ing dis- tance min	Focal length
	40.12	00	P	24.0	32.0
Achromatic	199.25	0.0	P	7.6	17.0
dry systems (45mm)	25/0.50	0	P	0.44	7.2
	400.65	D	P	0.26	4.6
Oil immersion objectives (45mm)	ÇI 1991.30	D	P	0.09	1.9
Flacrite	F1 46/0.85	01	P	3 0.20	3 42
dry systems 37mm)	F1 400.85	0	P	3 0.20	1 42
Fluorite oil immersion	FI (7) 54/0.96	DO	P	0.17	3.4
objective (37mm)	FI (7) 95/1.32	0.5	P	1 000	1 20
8, 9, 8	Iris FI CI 95/1.32 - 1.10	D	P	100	1 20
	Apo 12.5/0.30	DO	p	2.3	13.0
Apochromatic	Apo 25/0.85	D	P	0.76	7.4
dry systems (37mm)	Apo 40/0.95	019	P.	0.09	4.6
	Apo 63/0.95	019	P	0.09	3.0
Apochromatic oil immer-	Apo Cli 991.32	D	P	0.09	2.0
sion objectives (37mm)	Apo CI 991.40	D	P	0.03	2.0
	NP1 6.3/0.20	0.0	P	20	26.0
Plano objectives (45mm)	NP1 199.25	DO	P	0.53	16.0
NPI dor 18mm	NP1 19/0.40	D	P	0.50	11.0
field-of-view diamotor:	NPI 25/0.50	D	P	0.38	7.0
inter-co-sider distributed	NPI 495.85	0	P	0.15	4.5
	NPI OI 1001.30	D	. P.	0.26	1.7
	PI 1034				
	single-lens condensor	DO	P	30.0	33.0
Plano objectives (45mm)	PI 2.5/0.08	00		11.0	56.0
Plato cojectives (vomin)	PI FI 40.14 PI FI 100.30	00		14.0	40.0 18.0
Tor Money	P1 F1 150.30	00		1.26	16.0
(ield-of-view diameter)	PI 150.40 PI Apo 150.40	00	- 5	1.25	10.9
new-co-treat diaments	P1 Apo 19/0 40 P1 25/0.50			0.80	7.6
	P1 40 0.65		-	0.60	4.7
	Pl App 49 0.75		P	0.64	3.9
	P1 App GI 1001.32	0.5		0.74	2.4

glass thickness should be observed within ±0.05mm].

O: to be used without coverglass, DO: may be used with or without poverglass.

Di: coverglass thickness 0.17mm should be observed to an accuracy of +0.05mm, to set to this accuracy where the objective has a correction mount if it me.

9 Objectives in correction mount with automatic sharpness compensation. Its adjustment has hardly any effect on image sharpness, ideal method of focusing when the coverglass thickness is utknown, if These oil immension objectives may also be used for uncovered objects (sanear preparations without coverglass). The negligible reduction of image quality All I FITZ Plano, and Normal-Plano objectives have a total length of 45mm and are matched on the revolving nosepiece. This makes only negligible re-focusing with the fine adjustment necessary after a change of magnification.

The PI 1/0.04 objective with a total length of 65.6mm is an exception! This objective can be used only in association with the low-power condenser supplied with it and with the collector lens (1.13) in the foot of the stand swung-out of the optical path.

If special circumstances require the use of objectives of an adjustment length of 37mm together with those of 45mm adjustment length, their adjustment lengths must be increased to 45mm with the adapter Code No. 519164. The consequent increase in tube length is allowed for by a lens built into the adapter.

Oil immersion objectives are distinguished by their engraving . Oel" and a black ring around the bottom rim of the objective mount. It is a characteristic feature of immersion objectives that the refraction of the rays as they emerge

from the coverglass is reduced or completely eliminated and with larger angles of aperture the total reflection on the surface of the coverglass is also eliminated. This makes it possible for rays of larger angles of aperture to enter the microscope objective, which means an increase in the numerical aperture and therefore in the resolving power. The immersion oil has approximately the same refractive index n = 1.515 as the coverglass and the front lens of the microscope objective. Focal length and working distance of an immersion objective are usually very small. For this



Fig. 18 Complex Oel 100/1.30

reason care is indicated during work with such objectives. The coarse adjustment should be used only until the immersion objective has dipped into the oil: check by looking along the top of the coverglass. Focusing must be carried out with the fine adjustment only, and with continuous microscopic control. Ensure that the immersion oil is free from air bubbles. Use LEITZ immersion oil or, for fluorescence observation, non-fluorescent LEITZ immersion oil Generally the condenser No. 602 will be adequate also with oil immersion objectives, if, however, the full aperture of the immersion objective is to be used,

oil should also be introduced between the condenser top and the underside of the object slide. After the end of the investigation all obtical surfaces wetted with immersion oil must be carefully cleaned. For this purpose a soft rag spaked in xylene is suitable. Polish with a dry rag. Alcohol (methylated spirits) must never be used for cleaning the objectives and condensers. Avoid pressure during cleaning.

e.g. for very delicate structures, the

aplanatic-achromatic condenser No. 603.

A 1.25 should be used. Here immersion

In the optical system of the microscope. objective and eveniene form a unit. In the ORTHOLUX II field-of-view indices of up to 18mm can be achieved. The eveniene field of view is the area of the intermediate image in the tube that can

be surveyed with the eyepiece. It is magnified by the eveniene factor. The image diameter of a PERIPLAN GF 10 x eveniene of 18mm field-of-view diameter (field-of-view index 18) thus appears as large as a diameter of 10x18 = 180mm at a distance of 250mm from the observer. At this size the picture can also be reproduced on a groundglass screen 250mm above the eyepiece. If the diameter of the field-of-view is di-

vided by the objective magnification, the diameter of the object area that can be surveyed is obtained. With the above-mentioned PERIPLAN GF 10 x eyepiece and a 25/0.50 objec-

tive an object area of

= 0.75mm diam. can thus be surveyed.

Evenienes are generally constructed so that the observ's eye must be about 8 to 10mm above the top lens surface (evelens) (pupil distance). The so-called highpoint eyepieces have a pupil distance of about 20mm and therefore permit observation through spectacles and thus compensation of astigmatism. The highpoint eyepieces have a plastic top rim: this prevents the evenieces from being scratched.

The tables below offer a survey of the eyepieces that can be used in the ORTHOLUX II microscope:

DEDIDI AN eveniece

for scoepting a graticule.

Magnification	Field of view
6.3 x	18
6.3 x M	18
8 x	16
10 x °	14
25 x	8
High-point	10-10-318
8 x	18
10 x	15
10 x M	15
10 x MF1)	15

10mm = 100 intervals M - evepleces with focusing evelens and suitable MF - eveplece with focusing evelens and focusing Widefield eyepieces PERIPI AN NG/GE (diam. 23 2mm) Field of view Magnification

NF 10 x 18 NE 10 v M 18 GF 10 x 18 GF 10 x M GE 125 v GF 125 V M GF 125 x MF 1 GF 16 x GF 25 X 10 GE 25 v M 10

10mm = 100 intervals 5 with graticule for ORTHOMAT-W or for attachment camera with Pelaroid back or for system attachment

Special evenieces (diameter 23 2mm) Magnification Designation H 63x Pointer eveniece H 6x Pointer double evepiece P 10 v 0 Eyepiece with fixed pointer (° yellow engraving) NF 10 x * Eyepiece with fixed pointer (* yellow engraving) GF 10 x * Eveniece with fixed pointer (* vellow engraving) P 10 x Comparison eveniece GF 10 v * Evepiece for photomicrography (engraved red dot) 125 x Screw micrometer eyeplece 12.5 x Screw micrometer eveniege for monochromatic light Attach the microscope slide to the obtect stage.

Choose a medium-power objective, preferably NPI 10/0.25, with PERIPLAN GF 10 x eveniece for first examination.

Raise the swing-out condenser to its topmost position and swing the condenser top into the beam. Open the aperture and field diaphragms. Turn the swing-out lens in the foot of the stand into the beam by means of the knob: the lens remains in the beam during all examinations in brightfield and darkground - During work with the PI 1/0.04 objective this lens must be swung-out. Focus the specimen with coarse and fine

adjustment, Compensation for eyes of different visual acuity is carried out as follows:

Look through the eveniece with fixed evelens with one one eve (right hand evepiece/right eve), focus the specimen with the fine adjustment. Now look at the same area of the specimen with the other eye through the eyepiece with focusing evelens and adjust the latter until this area also appears sharp. -The fine adjustment must not be operated during this procedure. - After centration of the condenser, this adjustment must be exactly repeated and checked from time to time.

Now the field diaphragm must be centred (Fig. 19). The following rules apply to the use of

the field diaphragm and the aperture diaphragm:

The field diaphragm protects the specimen from unnecessary heat and prevents flare. It is therefore opened only far enough to fill the field of view of the

microscope objective.


The aperture diaphragm - as long as it is smaller than the aperture of the objective-determines resolving power and contrast of the microscopic image. Whereas with specimens of normal contrast range the aperture diaphragm is closed so that it only amounts to 2/3rds of the objective aperture, the following procedure is recommended with objects of low contrast range: the aperture diaphragm is opened far enough so that it is just visible in the rear lens of the objective (after removal of the eyepiece). The aperture of the condenser and that

of the objective are now idential. If at this setting all details of the specimen appear at sufficient clarity, the disphragm of the condenser is gradually closed until the less contrasty structural elements too become visible. It will in most cases be advisable to close the aperture diaphragm only far enough so that it transmits about 2/3rds of the full objective aperture. If it is closed further. the resolving power of the objective and with it the performance of the microscope rapidly deteriorate.

The aperture diaphragm must not be used for regulating image brightness. For this purpose only the transformer. or, with colour photography, neutral density filters should be used.

After the condenser has been centred, the specimen must now be fully illuminated. Remove the eyepiece from the eyepiece tube. Observing the rear focal

plane (visible aperture of the objective) adjust the setting knob of the lamp condenser until the entire specimen is evenly illuminated.

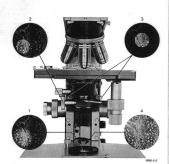

Centring the field diaphragm

Close the field disphragm
 By vertical adjustment of the condenser focus the image of the field disphragm visible in the

Move the image of the field disphraom into the

centre of the field of view with the two centring screws

Open the field diaphragm so that its rim just disappears become the field of view.

6 Transmitted-light-darkground

For investigations in transmitted-light darkground with the immersion darkground condenser D 1.20 mainly the special darkground objectives Oel 100/ 1.30-1.10 or FI Oal 95/1.32-1.10 with built-in iris diaphragm are used. In addition these immersion objectives can also be supplied without iris diaphragm: for darkground examination they must be used with a funnel stop to reduce their aperture, too high for the darkground condenser D 1.20, below the limiting aperture of the condenser (here 1.20) Otherwise part of the illuminating ray bundle would enter the objective and the darkground image would no

longer be perfect. Naturally, dry systems of large aperture can also be used with darkground condenser D 1.20. For darkground examination with dry systems of medium power, particularly

systems of medium power, particularly for serial examinations, the dry dark-ground condenser D 0.80, which is simpler to operate, is recommended. With dry systems of an aperture larger than about 0.70 an insert stop should be placed in the condenser 0.80. This stop, however, does not change the limiting aperture of the condenser; it merely absorbs stays light.

Adjusting the darkground image with the condenser D 1.20.

 Before inserting the darkground condenser set the centring mount roughly at the centre position (resetting with the two centring screwa). Push the darkground condenser into the dowald ground condenser into the dowald guide as far as it will go. Do not yet raise the condenser.

Place a drop of immersion oil, which must not be too small, on the condenser top surface.

 Place specimen on the object stage and focus it with the NPI 10/0.25 objective. If brightness is inadequate, lower

the condenser slightly, if necessary.

4) Raise the condenser, looking across its top, with the vertical adjustment until

the oil droplet touches the underside of the microscope slide (brief lighting up of the microscope slide). 5) Looking into the tube, move the con-

denser still closer to the microscope sidle, until the smallest possible sidle, until the sidle sid

6) Only now turn in an objective of higher primary magnification. Objectives of apertures larger than 1.15 should first be fitted with the appropriate funnel

7) When an immersion objective is used a drop of immersion oil must also be placed on the top of the coverglass. See also Use of oil immersion objectives in

transmitted-light-brightfield.

When the special darkground objectives are used maximum brightness is set with the fine adjustment of the microscope; at the same time the structure of the specimen must become weakly visible. Now close the disphragm on the objective until a perfect darkground image is produced: the object structures appear surrounded by a bright into

The various points concerning the adjustment of the darkground image logically apply also to the dry darkground condenser Do.B. Naturally immersion oil is used neither between condenser and microscope siden on between covergitas and objective. A suitable test object for darkground adjustment is a specimen of buccal spirilla, which can always be easily oblatined.

For details see Instructions No. 513-31.



On the right the centred light dot: the darkground condenser is correctly set



7 Investigations in phase contrast

For this purpose the phase contrast equipment according to Zernike is available. It consists of:

phase contrast condenser Phaco objectives Achr. + NPI auxiliary focusing magnifier two centring keys.

Operation of the phase contrast equipment:

equipment: Screw the Phaco objective into the revolving posepiece, and push this into the

changing guide; insert the phase contrast condenser, place the specimen on the object stage. Swing phase ring 1 in the phase contrast condenser into the beam. Swing in NP1 10/0.25 Phaco objective. Focus the specimen.

Focus the field diaphragm with condenser, and centre the condenser (see p.19 Fig.19). The aperture diaphragm remains fully open during all phase contrast investigations. Replace the evepieces with the facusing telescope. Adjust the eyelens of the telescope until the images of the light and the phase rings appear equally sharp. Check whether both rings are concentrically superimposed, if necessary establish coincidence by means of the centring keys. Replace the focusing telescope with the evepiece. If necessary, repeat the centring procedure after magnification change. For details see Instruction No. 517-84a.



Fig. 27 Phase ring and light ring as they appear in the focul ing magnifier, from left to right: a) in brightfield b) in phase contrast in centred position, c) in phas contrast in de-centred position, d) in derkorprovid




Phase contrast condenser in position 2 on ORTHOLUX II

Fluorescence

Phaco objectives

NPI Phaco objectives

Con- denser setting	Investigations in	Objectives
н	Brightfield	All objectives
1	Phase contrast	Phaco 190.25 NPI Phaco 190.25
1	Phase contrast	NP1 Phaco 160.40
2 2 2 2	Phase contrast	Phaco 250.50 Phaco 400.65 NPI Phaco 250.50 NPI Phaco 400.65
3	Phase contrast	Phaco 1001.30 NPI Phaco 1001.30
4	Darkground	All objectives

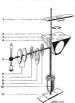

Suppression filter slide 9 EO suppression filter 3 Turnet with FO suppression filter and beam-splitting mirror, 4 Beam splitting mirror, 5 Light source. 6 Filter in the lamp housing, 7 Field disphragm, 8 Exciting-filter turnet, 9 Stray-light stop, 10 Objective. 11 Specimen

For the observation of specimens emitt-

the following outfits are available for the ORTHOLUX II microscope:

Transmitted-light fluorescence

ing primary and secondary fluorescence Mirror Housing 250 O with Lamp Housing 250 for 200W mercury lamp; dark-

ground condenser.

Incident-light fluorescence Mirror Housing 250 O with Lamo Hous-

ing 250 for 200W mercury lamp or Lamp Housing 100 Z with 100 W mercury lamp:

fluorescence vertical illuminator. The appropriate combinations for exciting and suppression filters are described

in the instructions for the Lamp Housing 250 and in the List Incident-light Fluorescence. Generally a dark background is aimed at in fluorescence microscopy The ideal filter combination has been found when these conditions are met. Directions: Switch on the lamp with the power unit and centre it according to instructions "Lamp Housing 250" No. 514-72b. Insert the exciting filter in the filter slot in the lamp housing, the suppression filter in the filter slot in the microscope.

Appropriate exciting filters in a filter turret are built into the fluorescence vertical illuminator Set the desired position Focus the specimen and observe fluorescence. If necessary change exciting and suppression filters.

Note: The brightness of the fluorescence image decreases as the square of the Eveniece magnification increases and increases with the aperture of the objective. Objectives of large aperture and evepieces of low magnification are there-

Detailed information about the fluorescence vertical illuminator can be ob-

fore recommended tained from the instructions No. 512-92

9 Investigations in incident-light darkground

With the ULTROPAK vertical illuminator the object is illuminated from all sides by means of ring condensers. Only by means of ring condensers. Only structure forms the lingsh diffusely reflected by the object structure forms the image. Observatoruler forms for indicated and carried out in incident-light dark countries of the images. Observatoruler forms described and object of the images. Observatoruler for the indicated on the images. Observatoruler for the indicated of the images of the images of the images of the images. Observatoruler for the elimination of disturbing reflections. Illumination of disturbing reflections. Illumination of irre-outlar and unoncernad surfaces.





Fig. 24 ULTROPAK vertical illuminato



The following ULTROPAK objectives are available:

Designation	of objectives Magnification/	Free working distance	*	ognificul th eyepi	ece	Outfits with immersion attachments on request
	aperture	in mm	63×	8×	10 x	0-110
	UO 4010	13.2	25 x	32 x	40 x	4
	UO 6.5/0.18	16.0	40 x	52 x	85 x	6.5
ry systems	UO 11/0:25	5.7	20 x	88 x	110 x	11
	UO 22/0.45	2.1	140 x	175 x	229 x	
	UO 33/8.55	1.0	200 x	255 X	330 x	
	UO 42/0.60	0.65	265 x	335 x	420 x	
mmersion	UO W 55/0.00	0.59	365 x	440 x	550 x	Dipping Cone EZ 23 - 100
bjectives	UO (01 60/0.85	0.36	380 x	480 x	600 x	together with 50 cover- glasses and 1 bottle of

Operation of the microscope for incident-light darkground investigation Attach lamp housing to the upper atta ment tube of the microscope according

Attach lamp housing to the upper attachment tube of the microscope according to instructions p. 5, Fig. 6. Insert the ULTROPAK into the horizontal

changing guide, Fig. 26, and tighten it with clamping screw (1.5). Insert the ULTROPAK objective.

Establish the best possible illumination by adjustment of the lamp condenser. With transparent specimens of for which the properties of the ULTROPAK where It can be ordated through 90°. The rear of the ULTROPAK commodates sector stops, for colour littles they additional softs will be properties of the properties of t

alyser is pushed into the filter slot (Fig. 1.4) of the stand.

Our instructions 513-90 give detailed information on further points, such as the use of the relief condenser, immersion attachments, and dipping cones.



Fig. 26 Inserting the ULTROPAK in the ORTHOLUX microscope

10 Measurements with the microscope

Linear measurements of specimens are normally carried out with a micrometer eyeplece* (graduation usually 10mm = 100 intervals), either during visual observation through the inclined tube of the binocular tube or on the focusing groundglass screen of the mirror reflex.

Before measurements are begun the micrometer value of the objective used must be known. The micrometer value is the length in the object plane which is precisely accommodated by the intervals of the graticule scale in the measuring eveniece after magnification by the objective. Although these micrometer values are laid down in tables, it is recommended to determine them with the aid of a stage micrometer, since the ontical constants of the objectives are subject to slight variations. The micrometer value remains the same also with photography with the large-format camera since the microscopic image and the graduation of the micrometer evepiece change at the same ratio as the bellows extension.



th image of the stage micrometer (on the right)

Example:

Determination of the micrometer value with the aid of a stage micrometer 2mm – 200 intervals and a micrometer open cown that gradual from — 100 intervals may be used to be used

12.20 μm. With low-power objectives, which do not form an image of the scale of the stage is of the stage in the stage of the stage is of the stage in the stage in the stage incometer across the stage incometer could be staged in the stage incometer coulded with 0 timerals in the microscope to stage and stage is staged in the stage incometer coulded with 0 timerals in the microscope that one of the stage incometer will be 0.38 in the 30 stage in the stage in the stage in the most accurate measurements under the microscope the screw microscope the sc

11 Heating stages

Heating and cooling stage 80

In place of the object stage No. 976, insert the rotating stage base No. 933. Turn the centring screws on the rotating stage fully back (without unscrewing them from the threaded hole). Push the object stage against the spring and lift it

out of the centring base.

Replace it with the heating stage. Further operation is described in our Instructions 515-81.

Heating stage 350 Heating stage 1350

These two heating stages can be used both with the object stage 976 and with the rotating stage No. 963.

Further details on the operation of the two heating stages should be obtained from the Instructions 515-89 and 515-74.



Heating stage 1350°

12 Photomicrography

General notes

These instructions have been written for the correct application and use of our ORTHOLUX II research microscope and cannot be regarded as a quide to photo-

micrography. We nevertheless want to point out a few important basic rules which must be followed without fail-

To obtain a good-quality photomicrograph, the following measures are

1) exact setting of Koehler's Illumination (para 4.3, p. 10), 2) the accurate focusing of the groundglass screen (large format) or in the eve-

piece or focusing telescope (35mm photography),

3) the accurate determination of the necessary exposure time,

4) meticulous cleanliness of all optical surfaces accessible to the user, such as deflecting mirror, dust glass, condenser, objective, tube lens, eveniece etc. (see section Care and Maintenance). Particular attention should be paid to

useful magnification, the choice of a light filter for the faithful reproduction (black-and-white photography) of the specimen.

the setting of the correct colour temperature of the low-voltage lamp for colour photography. and the choice of suitable exposure material for photomicrography.

Instruction No. 540-33

The following devices are available for photomicrography with the ORTHOLUX II-

ORTHOMAT® fully automatic microscope camera . . . Instruction No. 540-19 Micro-attachment for the I FICA® with vibration damper

System attachment camera In preparation CB 100 Polaroid* camera Instruction No. 540-34 9 x 12cm or 4 x 5in large-format camera] In combination with the 9 x 12cm or 4 x 5in large-format camera ARISTOPHOT® basic with fully automatic exposure control stand No 540_12 540_29

* Potaroid - Trademark of Polaroid GmbH

The range of applications of the ORTHO-LUX II can be extended with the accessories listed below:





Fig. 29 Discussion tube

Fig. 30 Comparison tube



Fig. 31 System attachment camera (with LEICA)



Fig. 32 Attachment camera with CB 100 Polaroid back

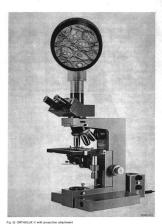


Fig. 33 ORTHOLUX II with projection attachment

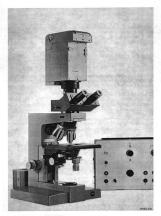


Fig. 34 ORTHOLUX II with ORTHOMAT W

13 Care and Maintenance

For protection the microscope is covered with the flexible dust cover after use. From time to time the stand should be cleaned with a linen or chamois-leather rag. Spirit must on no account be used. since it attacks the varnish. Petrol. however, is eminently suitable for the clan-

ing of varnished parts. Bright patches on the objective stage caused by petrol can be removed by

rubbing it with liquid paraffin or acidfree vaseline.

Special care is indicated during investigations involving acid (particularly acetic acid) or other corrosive chemicals. Direct contact of ontical surfaces and stands with these chemicals must be avoided at all costs and all parts should be thoroughly cleaned after use. The ontical parts of the microscope must be kept scrupulously clean. Dust on glass surfaces is removed with a fine. dry sable brush: this should be assisted by gently blowing across the glass surface at the same time. If the dirt resists removal, a piece of clean linen or a soft piece of chamois leather moistened slightly with distilled water should be used. If even this does not help, petrol or xviene, but under no circumstances alcohol (methylated spirit) should be used

For cleaning objectives should not be dismantled. If internal damage of objectives becomes evident these should be returned to the factory for renair

Special care is indicated during the cleaning of the reflex-reducing films. The external surfaces of the evenieces and the front lens of the objectives are coated with films about as hard plass. They should be cleaned with the same amount of care as uncoated glass surfaces. Some of the internal surfaces of the objectives and evenienes, however, are coated with very soft films, which must only be blown across with the greatest care. For this reason cleaning of the internal surfaces of evenieces is discouraged.

Careful maintenance preserves the performance of a LEITZ microscope for many years. If, however, an overhaul or repair of a damaged instrument is required our agencies or our factory should be contacted.

ensions and weld Height 414mm 253mm

Depth 253mm Weight 10,5 kg

14 Appendix

Magnification:

V= 25cm focal length objective or eyepiece in cm

Microscope magnification: V Microscope

V Microscope - M objective X V eyepiece

Magnification of the microscope with built-in tube lens:

V Microscope = M objective x V eyeplece x Tube factor

Photomicrography: Large-format camera

V Microscope bellows extension in cm

ORTHOMAT-W

M_{objective} x tube factor x eyepiece setting x camera factor 0.32

Projection tube:

At viewing distance 25cm = M objective X V eyepiece 15x

Projection prism

Vprojection magnification = V microscope x distance in m x 4

Visual projected image - projection magnification 4 x observation distance in m

Useful total magnification:

Vuseful magnification = 500 x A to 1000 x A

or

V_{eyepiece} = $\frac{500 \times A}{M_{objective}}$ to $\frac{1000 \times A}{M_{objective}}$

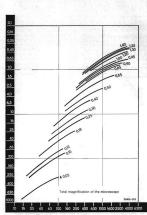
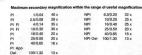


Fig. 35. Axial resolving power in μm for biological specimen n=1.45 (mean value) and $\lambda=550 nm$



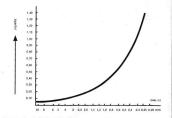



Fig. 36 Resolving power in λ and for green light (λ = 550nm = 0.55 μ m) as a function of the opertur

for the ORTHOLUX II microsc

Tube factor 1 x

Objective		1	otal magnifical	ion with eyepied	00	
magnification	6.3 x	8 x	10 x	12.5 x	16 x	25 x
1	6.3		10	12.5	16	25
2.5	16	20	25	32	40	63
3.5	22	28	35	45	55	90
4.	25	32	40	50	63	100
6.3	40	50	63	80	100	160
10	63	80	100	125	160	250
12.5	80	100	125	160	200	320
16	100	125	160	200	250	400
25	160	200	. 250	320	400	630
40	250	320	400	500	630	1000
54	340	430	540	675	860	1350
63	400	500	630	800	1000	1600
90	565	720	900	1125	1450	2250
95	600	760	950	1200	1525	2375
100	630	800	1000	1250	1600	2500
95	600	760	950	1200	1525	

ed-light beam in the ORTHOLUX II largefield microscope

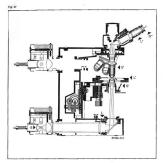

The beam puth is based on Koehler's principle of illumination. An image of the light source S is formed in the front focal plane of the substage condenser by the lamp collector iens. The aperture disphragm is also situated in this plane. The Image S' of the light source will then he reproduced by the condenser and the objective in the latter's rear focal

field diaphraom L is reproduced at a reduced scale in the object plane L' by the condenser. The objective intermediate image plane, and this image is once more viewed at a magnification through the eyepiece. more viewed at a magnification brough the eyepiece. The third image of the field diaphragm L'" appears together with the microscopic image on the retina plane S" and finally in the exit pupil of the eyepiece S". All these planes are optically conjugated. The



Registered Trademark
 Design subject to alteration without notice.

List 512 - 120 / Engl.

ERNST LEITZ GMBH D-6330 WETZLAR GERMANY Subsidiary: Ernst Leitz (Canada) Ltd., Midland, Ontario

WITH A VIE